

Notes for Analog Modulation

For the Analog modulation laser modules, please note that there are two lines (red+, black-) in the circuit panel for analog signal (0-5VDC) input.

1. Connect the laser head with the power supply correctly (Please note that the labels on the laser head and the power supply must be the same for ONE laser system. Unmatched connection will make the laser system fail.)
2. Without signal input (or the lines open), the laser is OFF.
3. With signal 0VDC input, the laser outputs Min value/No output.
4. With signal 5VDC input, the laser outputs Max value.
5. With other voltage between 0-5VDC, such as 1V, 2V, 3V, 4V, 4.5V, the laser outputs different powers.